



530 Main Street, Acton, MA 01720

Phone: (978)263-3584, **Fax:** (978)263-5086

Web Site: www.acton-research.com

**Supplemental Instructions
for
MF-460**

Ver 2.0

1
MF-460Supp.DOC

Per customer specification, this SpectraPro is supplied with an MF-460 slit adapter, which mounts to a SpectraPro entrance or exit slit, and accepts accessories with a 4.5” diameter bolt circle. To mount the MF-460 the following procedure is recommended:

1. Remove the Shipping Cover from the slit and attach the MF-460 to the slit using the provided four (4) #8-32 x 3/8 long socket head cap screws, assuring that the label “TOP” is on the top.
2. The MF-460 is now ready to accept an accessory with 1/40 –20 screws on a 4.5 diameter bolt circle.